
INDOOR
biotechnologies ^{INC}



DUSTREAM[®] Collector



Dust Collection and Extraction
in a Single Device

WWW.INBIO.COM



COLLECTING DUST

1. Insert the plastic filter into the DUSTREAM[®] collector and attach the collector to the vacuum cleaner tube.



DUSTREAM[®]
collector + filter



DUSTREAM[®] collector
+ adaptor



DUSTREAM[®] collector +

2. If the collector does not fit the vacuum cleaner tube, attach the adaptor piece to the collector.*
3. Turn on the vacuum cleaner and vacuum 4 separate areas of 8" x 12" (or approximate the size of a letter size paper) for 30 seconds each (total sampling time, 2 minutes; total area sampled ~0.25sq meter)



4. Remove the filter containing the dust sample and place it in a small Ziploc bag or envelope.



5. Rinse the DUSTREAM[®] collector with water, dry with a paper towel and place a clean filter inside for repeat sampling.

DUSTREAM[®] filter

6. Send the dust samples to Indoor Biotechnologies, Inc. for allergen analysis.

** Use the side of the adaptor which fits your vacuum cleaner.*



***DUSTREAM*[®] SAMPLE EXTRACTION**

- Two caps are provided with the *DUSTREAM*[®] collector system.
- Caps enable dust to be extracted with buffer within the collector and are designed for use with our rapid dust mite test.
- Caps can be used to prepare extracts on site if necessary, but **do not** need to be used for samples submitted to our *INDOOR*[®] Allergen Analysis Service.



DUSTREAM[®] Sample Extraction



SUBMITTING SAMPLES:

1. Fill out the INDOOR[®] Allergen Analysis Request Form completely and indicate which tests are requested.
2. Submit the samples to:

Indoor Biotechnologies, Inc.

INDOOR[®] Allergen Analysis Laboratory

700 Harris Street,

Charlottesville, VA 22903



DUSTREAM[®] Collector

- *Dust collection and extraction in a single device*
- *No dust handling necessary: Only 2 minutes sampling time*
- *Replaceable filters - test multiple sites with a single collector*
- *Universal adaptor—fits most vacuum cleaner models*
- *Easy seal leak proof caps*
- *For allergens, molds, endotoxins, other biologics*
- *Cost effective*



INDOOR® Allergen Analysis Request Form

Sample ID Sample Type Date Sampled Analysis Requested

Name

Address

Phone

E-mail

Method of Payment



Credit Card #

Exp. date

Signature

INDOOR

biotechnologies^{INC}



Indoor Biotechnologies, Inc.
700 Harris Street
Charlottesville, VA 22903

Phone: 434-984-2304

Fax: 434-984-2709

E-mail: mail@inbio.com